
MEMORANDUM




TO: DMT SUPERVISORS
FROM: AMANDA BOLDOC, ALCOHOL SECTION SUPERVISOR
SUBJECT: NEW DMT SIMULATOR CONFIGURATION
DATE: MARCH 23, 2017

Old Simulator Configuration:



The old simulator connected to the DMT via a tower with locking arms. The simulator head contained connections for the air flow and the temperature.

New Simulator Configuration:

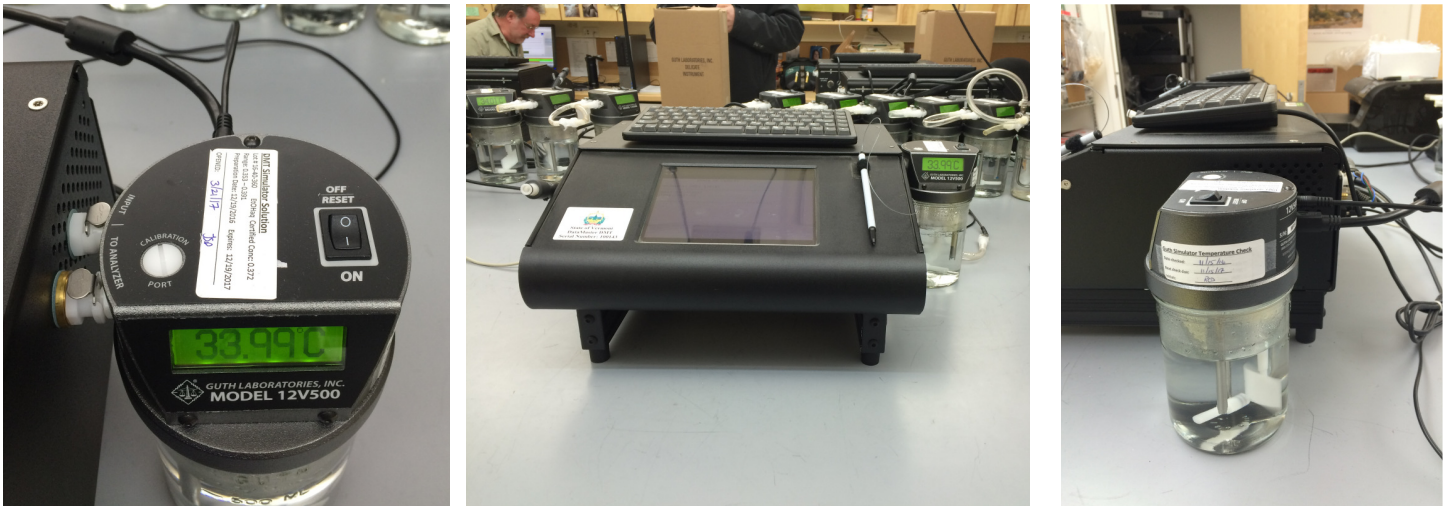

The new simulator directly connects to the side panel of the DMT via the air ports. The temperature cable connects to the DMT via an RS232 cable which is a small port on the simulator head and a large connection at the rear of the DMT. This simulator also has an on/off switch on the top. Software protocols (RPC and Simulator Solution Change) and maintenance procedures remain the same.

DO NOT PLACE ANYTHING UNDER THE DMT, NOT EVEN PAPER!

THE SIMULATOR MUST REMAIN ON AT ALL TIMES.

FOR QUESTIONS: amanda.bolduc@vermont.gov (802) 241-5300

Rob.driscoll@vermont.gov (802) 585-0123