

STATE OF VERMONT
DEPARTMENT OF PUBLIC SAFETY
VERMONT FORENSIC LABORATORY

P.O. BOX 47
WATERBURY, VERMONT 05676-0047




TEL: 802•244•8788
FAX: 802•241•5557
E-MAIL: cjs0lab@dps.state.vt.us
<http://www.dps.state.vt.us/cjs/>

MEMORANDUM

TO: DATAMASTER DMT SUPERVISORS; GREG NAGURNEY
FROM: TRISHA CONTI, SENIOR FORENSIC CHEMIST, ALCOHOL PROGRAM SUPERVISOR
SUBJECT: DMT HISTORY REPORTS
DATE: 8/9/2012

It has been observed that a “fatal error” may occur when **generating history reports** on the DataMaster DMTs. These errors will cause the instrument to reboot. This is due to a glitch in the software that is currently being used by all DataMaster DMTs in the State. We are aware of the situation and are working with the instrument manufacturer to rectify the problem.

When entering a date in the “End” date range box (after double tapping to highlight the field), you **MUST HIT “Backspace”** so the field shows MM/dd/yyyy. At this point you can enter in the desired date using the numeric keys.

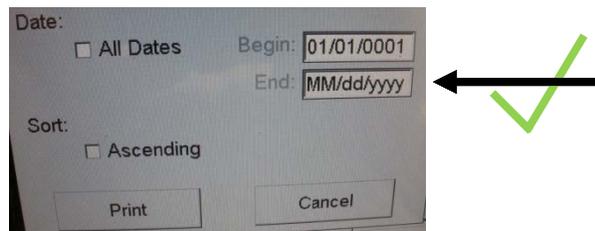

However, if you attempt to enter a date in the “End” date range box (after double tapping to highlight the field) without hitting “Backspace” first (so the field shows 12/31/9999), the DMT will encounter a “fatal error” and reboot.

